

CroPharm

Solutions for Pharmaceutical and Nutraceutical



Tube Coders

Tube Embossing Machines


www.cropharm.com

Coding Machines Type M & Type P

- Designed for single or double side coding of tubes with letter/ numbers on the fold of metal tubes or on the welded seam of plastic or laminated tubes. Rippling jaws for the seals are also available. With a built-in cutting device any irregularities of the welded seam of plastic or laminated tubes are neatly trimmed.
- Once the machine has been fitted with the required jaws it is adjusted for height according to tube length. Filled tubes in multiple tube holders are placed between the jaws of the machine.
- Coding and/or trimming is achieved by depressing a hand lever (COD manual) resp. by actuating the foot switch (COD Pneumatic). Correct positioning of each individual tube in the machine is ensured by a location peg which corresponds with holes in the tube holder.



COD M

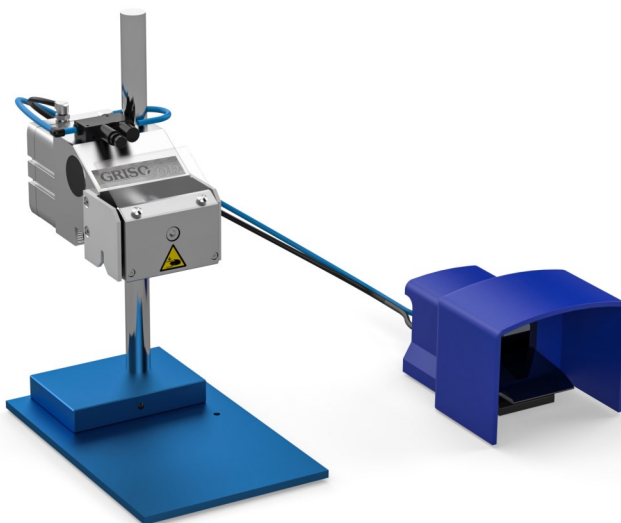


COD P

Technical Data		
Tube dimensions	length	20-320 mm max.
	Tube Ø	12-60 mm max.
Output	Up to 1500 tubes per hour	
Script Height	1.5 mm standard; (2 mm on request)	
Number of Digits	Max. 12 (1.5mm) 9 (2 mm)	
Working width	90 mm	
Dimensions	width	200mm
	depth	300mm
	height	420mm
Weight	approx.	8 kg manual, 11 kg pneumatic
Accessories	multiple tube holder, Set with coding digits	

Coding Machine Specifications

Jaws	Flat Jaw	Marking Jaw	Marking/ rippling Jaw	Rippling Jaw 1mm	Rippling Jaw 3mm	Cutting Device
Part Number	A.130010	A.130020	A.130024	A.130025	A.130030	A.130040
Trimming	P 2x					P
Coding Single Sided	P/M	P/M				
Coding Double Sided		M 2x (P)				
Coding/Rippling	P	P				P
Coding/rippling/ crimping			P (set)			P
Coding/rippling			M (set)			
Crimping				P (pair)	M (Pair)	
Crimping Rippling				P (Pair)		P

Plastic (P) Metal (M)

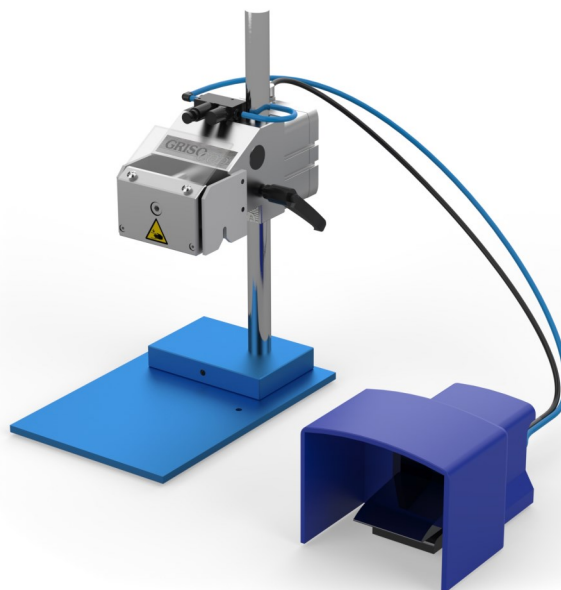



About Us

CroPharm is the successor of Chemical and Pharmaceutical Industry Co. Ramsey, NJ founded in 1956. We specialize in tablet dedusting and capsule polishing for pharmaceutical processing and production. CroPharm also offers solutions for tube filling, crimping, capsule filling, and tooling inspection. We offer engineering service, customization, and after sales support.



@cropharm.inc



@CroPharm Inc



@CroPharm

Contents within this brochure are subject to change without notice.

CroPharm

132 Pepes Farms Rd
Milford, CT 06460

Call: +1 (203) 877-3859

Email: sales@cropharm.com

www.cropharm.com